Haier



Ultra Low Temperature(ULT)Freezer Operation Manual



Manufacturer: Qingdao Haier Biomedical Co., Ltd. Address: Haier Industrial Park, Economic Technology Development Zone, Qingdao P.R. China Web:www.haiermedical.com Revision Date:08/2022 Version:3rd,2022 Dedicated code:0270501129A V13026



Model DW-86L419 DW-86L579 DW-86L729 DW-86L829

- Read the Operation Manual carefully before using your appliance.
- Keep the Operation Manual in a safe place.
- Appearance , color and layout of the door may vary.
- Translation of the original instruction.

Content

Safety Labels and Safety Precautions	1
Usage Precautions	5
Product Feature	7
Installation	8
Freezer parts and control panels	12
Application method	14
Alarm	30
Cleaning maintenance	31
Recycling the Rechargeable Battery	33
Optional Accessories	34
FAQ	36
Refrigeration and Wiring Diagram	37
Specifications, Technical data and Packing list	39

Technical Data

Model	Net Volume (L)	Rated Voltage (VAC)	Rated Frequency (Hz)	Rated Power (W)	Weight (kg)	Dimensions (W×D×H) (mm)
DW-86L419	419	220-240~	50	800	255	830x980x1980
DW-86L579	579	220-240~	50	1000	300	903x980x1960
DW-86L729	729	220-240~	50	1700	350	1049x980x1980
DW-86L829	829	220-240~	50	1900	380	1153x980x1980

Packing List

Name	User Manual	Instruction to install sapcers	Plastic bag	lce Scraper	Key	Spacer
DW-86L419	1	1	1	1	4	2
DW-86L579	1	1	1	1	4	2
DW-86L729	1	1	1	1	4	2
DW-86L829	1	1	1	1	4	2

Global Warming Potential

Model	Rated voltage(VAC)	Rated frequency (Hz)	CO ₂ equivalent(Tonnes)
DW-86L419	220-240~	50	0.0036
DW-86L579	220-240~	50	0.0041
DW-86L729	220-240~	50	0.0048
DW-86L829	220-240~	50	0.0051

This product contains fluorinated greenhouse gases covered by the Kyoto Protocol. Do not vent into the atmosphere.

GWP=global warming potential

Refrigerant type	GWP
R170	20
R290	20

Specifications, Technical data and Packing list

Specification


Name	Ultra Low Temperature Freezer
Model	DW-86L419/579/729/829
Exterior/Interior wall material	Coated cold rolled steel
Outer doors	Coated cold rolled steel
Inner doors	Plastic framed PS board
Shelves	Stainless shelves(height adjustable)
Porthole for testing	2
Insulation	Vacuum insulated with polyurethane foam(non-CFC)
Compressors	High stage:hermatically sealed Loe stage:hermatically sealed
Evaporator	Copper tube
Condensor	Finned coil
Refrigerant	R170 R290
Temperature controller	Microprocessor controller
Temperature display	LCD display
Temperature sensor	RTD(Pt100)
Alarm device	High/Low temperature alarm, probe failure alarm, Hot condenser alarm, Ambient temperature alarm, Low battery alarm, Door ajar alarm, Power failure alarm
Remote alarm terminal & Battery	Maximum load:30V DC, 2A Rechargeable battery:12V DC, charges automatically DW-86L419/579/729/829:Lead-acid battery
Electric shock protection type	I
Noise	47dB(A)/48dB(A)/52dB(A)/48dB(A)
Ambinet temperature	10 to 32
Freezer temperature	-40 to -86
Foaming cabinet	HFO-1233zde
USB	Standard
RS485	Standard
pollution degree	2
overvoltage category	Ш
	40

Safety Labels and Safety Precautions

Dear Haier Customers,

Thank you for purchasing a Haier ULT freezer. Please read this manual carefully to avoid injuries to personnel and damage to the product. Please also follow descriptions marked with the following labels.




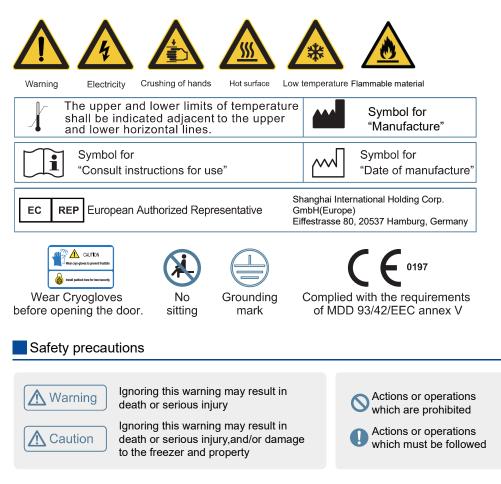

- When a CO₂/LN₂ backup system is activated, the installation place must be well ventilated. Increased CO₂ in the air may be harmful and even fatal. If the ventilation is poor, other methods should be considered in order to ensure safe working environments.
- If there is a leakage of petroleum gas or other flammable gas, close the gas supply valve and open doors and windows to ventilate the air. Do not plug or unplug your freezer unit in order to avoid potential explosion or fire.
- Only professional technicians or Haier service personnel can install the unit. Failure to do so may cause electricity or fire.
- The freezer must be securely installed on a firm floor. Tilted installation may result in the product tipping over thereby causing injury and damage.
- Please use a dedicated power supply matching the requirement specified on the label to avoid fire and shock.
- If the voltage being used is 10% higher than the rated voltage, a regulator with a capacity of 4000 W or higher must be installed.
- If the power cord needs to be extended, the cross-section of the extended cable must be no less than 2 mm² and no longer than 3 m for products of 220V~240V/50Hz to avoid fire or electric shock.
- Your Haier ULT unit is equipped with a standard three-prong power plug(grounded) complying with the standard three-prong socket (grounded) rated 16 A (220V~240V/50Hz). Removal of the ground prong is strictly prohibited for safety reasons under any circumstances. The electrical power plug should be securely plugged into the socket. A loose plug in the socket may cause fire.
- The power socket intended for your Haier ULT usage must be grounded to avoid electric shock.
- If the socket does not meet this requirement, the condition must be corrected by a qualified technician before using the ULT unit.
- Replacement of any spare parts should be performed by qualified technicians.
- Never install your ULT in an unprotected area. If the unit is rained on, there is a danger of electric shock.
- Your Haier ULT must not be installed in a damp area or an area subjected to water spray. Otherwise this may reduce the degree of insulation and thereby cause electrical leakage or electrical shock.
- 🚫 Never directly pour water into the unit. The water may cause electric shock or short circuit.
- O not place any water container or heavy object on top of the unit. A falling object may injure an operator. If the water spills into the unit, it may damage the insulation thereby causing electric shock.
- Never use gas lines, water mains, telephone lines or lightening rods as the grounding device for your Haier ULT unit. This type of improper grounding may cause electric shock or other danger.
- O not touch any electrical components, switches or power cord with wet hands. Such action may lead to electric shock.


Wiring diagram

•DW-86L419/579/729/829

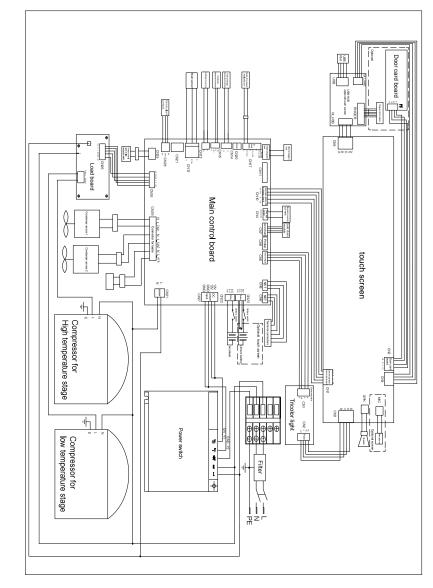

Refrigeration and Wiring Diagram

Refrigeration diagram

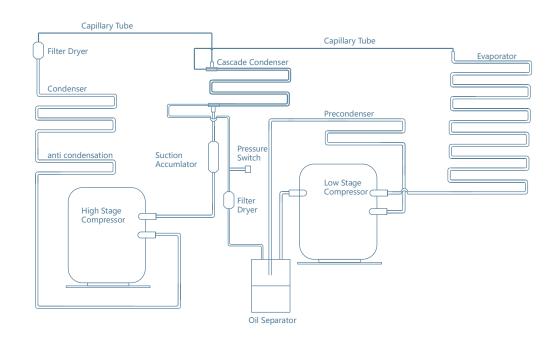

- When unplugging the power cord from the socket, please grip the plug itself and pull it out. Do not pull the power cord as this may strip the wires out of the plug thereby causing electric shock and fire.
- Should there be any malfunction in the equipment, power off the unit and unplug the power cord from the power supply. Continuous operation in an abnormal condition may result in electric shock and fire.
- Never disassemble , repair , or modify the unit yourself . Any such work carried out by an unauthorized person may result in fire , or electric shock or injury due to a malfunction .
- **D** Before any repair and maintenance of the freezer, please disconnect the power to avoid electric shock or injury to personnel.
- When repairing and maintaining your freezer, take precautions not to inhale any chemicals or aerosols floating inside and outside the unit. They might be harmful to your health.
- If poisonous, radioactive or other harmful materials need to be stored in the unit, the equipment should be located in a safe zone. Improper usage of the equipment with such materials may harm the environment or operator's health.
- If the unit is not in use for a long period of time, make sure the power cord is unplugged. Deteriorated insulation of the power cord may lead to electric shock or fire.
- If the unit is to be stored unused in an unsupervised area for an extended period , ensure that children do not have access and that doors are locked completely with a key.
- **D** The disposal of the unit should be accomplished by appropriate personnel . Remove doors to prevent accidents such as suffocation .
- 🚫 Do not use any non manufacturer-approved electrical components in the freezer.
- Never store flammable, explosive or volatile materials in the unit and do not use any flammable spray near the unit, as this may cause an explosion or fire.
- Never store corrosive chemicals with acid or alkaline properties in the unit as this can lead to damage to internal components of the unit.
- O Do not place any glass container or enclosed container into the freezer. These containers may crack at low temperatures causing injury to operators
- 🚫 Do not put the packing plastic bag within reach of children as suffocation may result .
- O not climb on top of the unit or place any object on it. Falling equipment may cause injury or property damage.
- O Do not use any hard objects such as nails and wires to explore any openings or gaps such as air ventilation ports. Accidental contact between a hard object and a moving part may result in electric shock or injury.
- Do not use electrical appliances inside the chamber of the appliance unless they are of the type recommended by the manufacturer .
- The appliance must be positioned so that the plug is accessible.
- The power plug is used as the disconnecting device.
- The appliance must be placed on a solid and flat surface, or excessive vibration and noise may be produced when the appliance in operation.

FAQ



Should there be any malfunctions in the system, please attempt to answer the following questions before notifying maintenance or calling a Haier Equipment & Instrument Service Center. Please do not dismantle the freezer yourself.

Fault	Troubleshooting
	Is the power supply normal? Has the main protector been turned on yet?
Freezer does not start up	Is the voltage supply too low?
	Is there any voltage input from the outside?
	Is the ambient temperature too high?
	Are the inner doors and outer doors closed properly? (Has any ice or frost damaged the seals between the door and the frame?)
	Is the condenser filter clogged?
Poor	Is the temperature setting correct?
refrigeration effect	Is the freezer being kept away from direct exposure to sunlight?
	Is the freezer near any heat source?
	Is the porthole plug installed in the porthole with proper insulation materials?
	Has the freezer been loaded with too many non-frozen samples within the last few hours?
	Is the unit set on a firm and level floor?
The unit is noisy	Is the exterior of the unit touching any objects?
-	Is the freezer unit leveled with the leveling legs?

- The applicance can be used by the persons with reduced physical sensory or mental capabilities or lack of experience and knowledge only if they have been given supervision or instruction concerning use of the appliance in a safe way and understand the hazards involved.
- If the power cord is damaged, it must be replaced by a qualified technician to avoid an electrical hazard.
- CP/IP foaming materials are flammable. They should be handled professionally.
- To avoid the risk of electric shock, this equipment must only be connected to a supply main with protective earth.
- There should be at least 30 cm space between the surrounding walls and the freezer for ventilation.
- Equipment cannot run in the condition of rich O₂ and flammable gas or liquid.
- After restarting your unit from a power outage or shutdown, ensure that all settings are correct. Accidental changes in settings may damage the stored products.
- In the event of a power outage and recovery, be sure to wait for at least 5 minutes before turning the unit on again to avoid damage to the compressors and refrigeration system.
- The air filter for the condenser should be cleaned regularly. A dirty filter could cause a malfunction or the freezer temperature to rise.
- During any repair operations, gloves should be worn to prevent getting injured by sharp edges or corners.
- O Do not use bare hands to directly handle any stored products. The cold temperature of the products and the interior walls may cause frostbite.
- Hold firmly onto the handle to close the door to avoid pinching your hands.
- O not tilt the unit more than 45 degrees when moving the unit.
- When moving the unit, please be careful not to stumble with the unit which could cause injury to personnel and damage to the unit.
- O not attempt to use the handle to lift or move the unit to avoid damaging the freezer or injuring personnel.
- Please open the lock first, then lift the handle.
- Maximum loading on the each shelf should be no more than 50 kg and total loading for whole unit should be no more than 200 kg. Heavier loads may cause damage to the shelving system.
- Keep ventilation openings, in the appliance enclosure or in the built-in structure, dear of obstructions.
- O Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended but the manufacturer.
- O not damage the refrigerant circuit.
- Unauthorized opening of the top cover of the control cabinet is prohibited in order to prevent damage to the inside components or injury to the operator.
- Turn the battery switch on before starting the unit, do not arbitrarily turn it off.
- When the ULT unit has been placed in a storage or not in use for a long time, its battery should be tested for low capacity because the battery may have already released all of its energy. Should this occur, please turn on the battery switch and run the unit for about a week to fully charge up the battery.

• If the freezer operates in authorized mode, first ensure that the magnetic card or fingerprint is registered on it. When you can create a new account, the registration will be completed by swiping the card or punching the fingerprint, meanwhile a magnetic card shall be assigned to the new account. Only when the magnetic card corresponds to the login account, can the electromagnetic lock be unlocked.




• If the user does not use the electromagnetic lock temporarily, it can be fixed with a clip, so as to ensure that the door can be opened / closed without swiping the card or using the fingerprint.

When there is a loss of power or the backup battery power is out of power, please contact technical support personnel should there be aa need to open the freezer door.
If the card is not swiped or fingerprint information is not entered, the door should not be opened forcibly as it can damage the electromagnetic lock.

5. Frozen storage rack and frozen storage box

When the freezer stores small size of materials, the frozen storage rack and the frozen storage box can be used, so as to utilize the space more effectively.

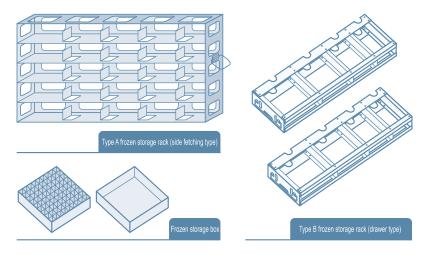

Model	Storage Rack(Type	Storage Box	
Widdei	variety	Amount	Amount
DW-86L419	5×5	12	300
DW-86L579	5×5	16	400
DW-86L729	5×5	20	500
DW-86L829	5×5	24	600

- Keep all ventilation openings in the enclosure or, in the structure for building-in, clear of obstruction.
- Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.
- Do not damage the refrigerant circuit.
- In order to reduce flammability HAZARDS the installation of this equipment shall only be carried out by a suitably qualified person.

Usage Precautions

1.When the unit operates normally, the door frame is slightly warm. The embedded hot gas halo tube provides free heat to prevent condensation to from on the door frame.

2.Before samples are loaded into a freezer, make sure the unit has reached the set temperature. Samples should be loaded into the freezer in batches. Each batch should be no more than 1/3 of the freezer capacity. This process allows the freezer to pulldown the temperature in a reasonable time.

3. The temperature display indicates the temperature where the temperature sensor is mounted inside the unit chamber, which may vary from the temperature at the center of the freezer, but it will gradually reach the actual temperature of the freezer over time.

4.Two access ports are available for testing in the back of the freezer. Thermocouple wires can go through the port holes to reach the interior of the freezer for temperature mapping. The gap of the holes should be sealed with insulation materials so that the interior temperature is not influenced by the ambient temperature.

5.When cleaning the unit, mild and neutral detergent solution should be used. Never use a hard wire brush, acid, gasoline, detergent powder, polished powder or hot water for cleaning. These tools and materials can damage the paint and coating of the unit. Particularly, never use gasoline or a solution with volatile chemicals to clean plastic or rubber parts.

6.After the freezer runs for some time, a layer of frost usually forms on the interior wall and inner doors. When the frost gets too thick, the refrigeration effect can be impaired. Energy consumption will increase. If the thickness reaches 5 mm, please use the supplied scrapper to remove the frost.

7.Before removing the frost, temporarily transfer the stored samples to another freezer. This is so that the temperature does not rise in the unit and damage the samples.

8.Behind the interior walls, there are many refrigeration tubes. Do not use a knife, an ice pick, or a screwdriver to cut ice and frost. This may damage not only the liner but also the refrigeration tubes.

9.If the freezer is not in use for a long time, please turn off the power and switch off the backup battery. The power cord should be unplugged.

10.Federal Communications Commission (FCC) Interference Statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules.

These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generate, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications.

However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one of the following measures:

•Reorient or relocate the receiving antenna.

•Increase the separation between the equipment and receiver.

Optional Accessories

1.Temperature recorder

When using the temperature recorder, please refer to the "User manual for Temperature Recorder" provided with the recorder.



tution The temperature recorder should only be installed by professionals or Haier serviceman.



ng Before installing the temperature recorder, please cut off the power supply to avoid electric shock or fire.

2.LCO₂ backup cooling system

For installation and operation of the backup cooling system, please refer to the user's manual provided with the system.



For LCO_2 back up system, please make sure the LCO_2 supply bottle has an internal liquid dip tube to make sure the liquid is fed to the backup system. For LN_2 backup system, please use a LN_2 dewar of 35 to 50 psig.



• Whenever a CO₂ or LN₂ backup system is installed, the location of the freezer must be well ventilated. Increased concentration of CO₂ in the air is harmful and even fatal. If the ventilation is poor, alternative methods should be considered to reduce the nitrogen or CO₂ concentration to the normal level.

• If a LCO₂ or a LN₂ cylinder fell to the ground, its valves nay get damaged, resulting in a risky condition. Fastening the bottles with a chain are to secure the them to ensure the bottles are positioned vertically and securely.

- The temperature of liquid CO_2/LN_2 is extremely low, which could cause frostbite. When replacing the cylinder, please always wear a pair pf protective glasses and protective clothes.
- This kind of backup cooling system should be used with the LCD panel.

3.Wireless cold chain module

The wireless cold chain module adopts wireless communication as a medium to realize the exchange between the specimen bank inside ultra-low temperature freezer and the specimen bank system, as well as the functions of data synchronization, management synchronization, operation synchronization and information synchronization.



4.Electromagnetic lock

• The electromagnetic lock can effectively enhance the safety of specimen access inside the freezer, which is convenient for use. When using it, make sure to swipe the card or use the fingerprint to unlock the lock before opening the door. If not swipe the card or use the fingerprint, or pull the handle forcibly, it may destroy the electromagnetic lock.

• Usage method: Before using the electromagnetic lock for the first time after unpacking, firstly open the door, and then remove the small clip for fixing the spring of electromagnetic lock.

• When the electromagnetic lock is unlocked, it may continuously maintain the open state for about 10 seconds. If more than 10 seconds later, you need to open the door once again, please swipe the card or punch the fingerprint. The unlocked/ locked state of electromagnetic lock is displayed on the touch screen.

• If the freezer operates in normal mode, any attached magnetic card can be used to open the door by swiping it in the card reading area of the display panel cover,but the fingerprint does not have the function.

Recycling the Rechargeable Battery

The Haier freezer is equipped with a rechargeable battery. This battery is recyclable. When the battery reaches the end of its life, please contact a local recycling organization for inspection or properly discard the battery.


1.Location of the battery

The battery in the electric cabinet is for the power outage alarm. It is located inside the control box of the right side of the unit.

To prevent electrical shock, only a qualified technician should replace the battery.

2.Removal of the battery

- •Turn off the power of the unit and unplug the power cord from the socket.
- •Use a screwdriver to remove the screws on the side panel and take down the side panel.
- •Unplug the connecting terminals from the battery.
- •Remove the bracket that fastens the battery Remove the battery.
- •Follow regulations to recycle the battery or discard it properly.
- When changing the battery, you must make sure that the brown wire connects to the positive pole of the battery, and the blue wire connects to the negative pole of the battery. The polarity must not be reversed. Incorrect polarity can damage the main control board so that it cannot charge the battery.

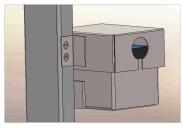

▲ Caution

After interruption of power supply, the battery can maintain the standby time of the touch screen for 22 hours.

•Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.

•Consult the dealer or an experienced radio/TV technician for help.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

(1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

FCC Caution: Any changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate this equipment.

11.RF exposure warning

This equipment complies with FCC radiation exposure limits set forth for an uncontrolled environment. This equipment must be installed and operated in accordance with provided instructions and the antenna(s) used for this transmitter must be installed to provide a separation distance of at least 20 cm from all persons and must not be collocated or operating in conjunction with any other antenna or transmitter.

This equipment complies with ISED RSS-102 radiation exposure limits set forth for an uncontrolled environment. This equipment should be installed and operated with a minimum distance of 20cm between the radiator and any part of your body.

Pour se conformer aux exigences de conformité CNR 102 RF exposition, une distance de séparation d'au moins 20 cm doit être maintenue entre l'antenne de cet appareil ettoutes les personnes.

12.Contains FCC ID: 2ASR8NFCV3, 2ASR8SC3832

13. This device contains licence-exempt transmitter(s)/receiver(s) that comply with Innovation, Science and Economic Development Canada's licence-exempt RSS(s). Operation is subject to the following two conditions:

(1) This device may not cause interference.

(2) This device must accept any interference, including interference that may cause undesired operation of the device.

L'émetteur/récepteur exempt de licence contenu dans le présent appareil est conforme aux CNR d'Innovation, Sciences et Développement économique Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes :

(1) L'appareil ne doit pas produire de brouillage;

(2) L'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.

Meaning of crossed –out wheeled dustbin:

X

Do not dispose of electrical appliances as unsorted municipal waste, use separate collection facilities. Contact you local government for information regarding the collection systems available. If electrical appliances are disposed of in landfills or dumps, hazardous substances can leak into the groundwater and get into the food chain, damaging your health and well-being.

Product Feature

This product is suitable for the ultra low temperature storage of products in applications such as clinical, pharmaceutical, scientific research, and epidemic institutions. It also can be used in blood stations, hospitals, centers for disease prevention and control, science and research institutions, electronic and chemical laboratories, biomedical engineering institutions, and open sea fishery companies to store red blood cells, viruses, germs, skin, bones, bacteria, sperm, biological products, electronic components, and low temperature testing samples of special products, etc.

1.Temperature control

Temperature is controlled by computer and numerically displayed, and regulated in units of 1°C (LCD:0.1°C); temperature range: -40°C to -86°C .

2.Security system

•Various malfunction alarms (high/low temperature alarm, power failure alarm, probe failure alarm,hot condenser alarm, high ambient temperature alarm, door ajar alarm, low battery alarm).

•Two types of alarms (buzzer sounding alarm, flashing light alarm).

•Multiple levels of protection are standard including passcode and time-delayed start.

•All components are electrically grounded.

3.Refrigeration system

•Optimized multiple refrigeration technology with top brand compressors off ering better refrigeration capability.

•Excellent temperature preservation layer providing optimal temperature preservation effect.

•Exclusive sealing structure of multiple doors and hot tubing for condensation prevention can reduce the frost effectively.

•Specially designed low temperature computer control, to prevent the normal redundant systems from being erroneously controlled by the low temperature compressor.

4.User-friendly Design

•Display to indicate freezer temperature, ambient temperature and input voltage. Control panel to set the high/low temperature alarm and inner temperature. All malfunction alarms are also shown on the display.

•Designed with adjustable shelves, suitable for product storage.

•Safe lock design prevents accidental opening of doors.

•Installation in broad range of ambient temperature range of 10℃ to 32℃ environment.

•Unique all-in-one latch design and compact caster features allow ease of operation and maneuvering.

•Condensing fan motors are automatically controlled to turn on and off for energy saving. •Network and remote alarm contacts are available as well for convenient connection and communication.

Due to our continuous improvement, Haier ULT freezers may appear to be different over time. User manuals are subject to change without notice.

·Load the samples back into the unit after it reaches the set temperature.

•Turn on the backup refrigeration system if there is one.



▲ Caution | Do not use any sharp tools such as knives or screwdrivers to defrost.

4.Battery maintenance

• When the control panel shows an alarm signal for "Low Battery ", please make sure the battery switch is turned on, and the battery will be charged. After about one week of charging, please recheck the battery capacity. If normal, the battery should be in full capacity. However, if the capacity is still low. please change the battery.

• The battery that supports the power outage alarm is a consumable item. The life expectancy for the lithium battery is 5 years(Lead-acid battery is 2-3 years.). If the battery is more than 5 years old(Lead-acid battery is 2-3 years.), the battery should be replaced because the alarm function may not work properly. To do so, please contact an Haier Equipment and Instrument Service Center.

5.Disposal of the freezer



• If the unit is to be stored unused in an unsupervised area for an extended period, ensure that children do not have access and that doors are locked completely with a key.

• The disposal of the unit should be accomplished by appropriate personnel . Remove doors to prevent accidents such as suffocation .

Cleaning maintenance



• To prevent electric shock or injury to operators, the AC power supply to the freezer must be disconnected completely before any repair and maintenance work is to be performed.

• During any repair maintenance work, do not inhale medical particles or aerosols near the equipment as they might be harmful to your health.

1.Cleaning the freezer

· Clean the unit once a month. This can help the exterior look new.

• Use a dry cloth to wipe away loose dust inside and outside of the freezer. If the unit is rather dirty, use a clean doth soaked with a neutral detergent to clean the unit. Then use a dry cloth to wipe away any residual detergent solution.

• Never pour water onto or into the unit. Doing so can damage the electric insulation and cause failure.

• Compressors and other mechanical parts are hermetically sealed. They do not need lubrication. Users should remove the frost on the chamber and clean the condenser filter as often as possible.

2.Cleaning the condenser filter

Clean the condensor filter when the control panel shows an signal for "Hot Condensor" and the alarm flashes. Even if the light is not on, the condensor filter should be checked regularly according to the suggestion from the distributor.

- To clean the filter, follow the procedure below.
- 1. Pull off the front grill cover.
- 2. Pull out the filter screen.
- 3. Use water to wash the filter screen.

4. After the filter screen is dry, reinstall it back in its original position and close the cover.

5. If the "Hot Condenser" light is on before cleaning, check the light to make sure that it shuts off after cleaning. If it does not shut off, please contact after-sales service personnel.

3.Defrost

Frost and ice can form in between the door gasket and frame to form an air gap, which can decrease the refrigeration effect of the unit. Please use the provided plastic scraper to defrost the interior doors.

To defrost the units, please refer to the instruction as follows.


•Turn off any backup refrigeration system if there is one.


•Remove the samples from the unit that needs to be defrosted. Move them to another unit or a container for temporary storage.

Turn off the power supply.

•Open the outer door and inner doors to let the unit thaw for a period of time.

•Use a dry doth to soak up and remove any water on the floor of the unit.

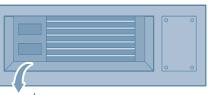

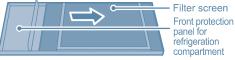

Installation

Installation environment

•Ambient temperature:10 $^{\circ}$ to 32 $^{\circ}$.The ideal temperature is 18 $^{\circ}$ to 25 $^{\circ}$ C. If necessary, use an airconditioning system to achieve the required ambient condition.

Ambient humidity: less than 80%. At an environment of 32°C, humidity should be less than 57% RH.
The intended location should be low in dust count.

- •The intended location should be vibration and shock free.
- •The highest elevation the unit can work safely: 2.000 m above sea level.
- •Input voltage: within Rated Voltage ±10 % .
- Overvoltage category: The transient state is Category II facility.
- Pollution Degree:2.



•An ULT freezer is usually sensitive to its operating environment. If a unit is not

installed in the conditions mentioned above, it cannot operate reliably. Please improve the environmental conditions before using the equipment.

•It is prohibited to use the unit in an outdoor place. After the unit is rained on, there is a danger of electric shock.

Installation site

For the equipment to achieve optimal operating conditions, an intended installation location should satisfy the following requirements.

•Do not install the unit in a confined place. The doorway should be large enough for the unit to freely enter or exit the room if necessary. This is to allow the unit to be repaired easily and timely to avoid damage to property

•The location for installation should be flat and firm.

•There should be good ventilation and no direct sunlight.

•The freezer unit cannot share the same power socket with other equipment. The power plug should be securely connected with the power socket.

•The power cord for the freezer should not be twisted or pinched.

•If the power cord needs to be extended, the cross-section of the extended cable must be no less than 2 mm² and no longer than 3 m.

•Before using the freezer, check the voltage supply. A voltage stabilizer to deliver rated voltage ±10% is recommended for areas where the voltage is known to be unstable. The voltage stabilizer should be rated at least 4000W.

•The freezer must be securely grounded

•If the power socket is connected with a ground terminal, make sure to inspect it for proper connection before using the equipment.

•If the socket is not grounded, it must be connected to a grounded terminal by a qualified technician.



•Do not ground the freezer through gas lines, water mains, telephone lines and lightning rods as this may lead to electric shock.

•After installation, the power plug must be easy to reach. In case of emergency, it is easy to unplug. Nothing should block the ventilation port of the freezer.

Preparation before Use

1. Remove all packaging materials, including base

Remove all the packaging materials, including base, at the right place, and remove the fixed tape or other attachments on the box for safe transportation.



A fixed connection strap is used between the left and right sides of the freezer and the base. Please remove the connection strap before removing the base;

• To handle the freezer before unpackaging, forklift or package carrier may be used. If the forklift is used, it shall be inserted at the bottom of blocking from the front or back of the freezer to move the freezer. If the package carrier is used, it shall pick up at the bottom of blocking and can only pick up the side of the freezer.

- After unpackaging, to handle the freezer, you may push wheels for handling.
- During handling, the inclination angle of freezer shall not be greater than 45°.
- 2. Check the supplied accessories

Check the items in the packing box according to the packing list. If they do not match each other, please contact Haier immediately.


3. Installation environment

There should be at least 30 cm space between the surrounding walls and the freezer for ventilation.

4. Adjust support legs

Rotate the leveling legs clockwise to extend them out to support the unit to the floor to ensure that the unit does not move while in usage.



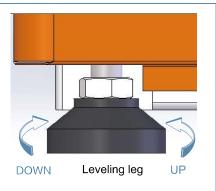

5. Placement

After adjusting and cleaning the unit, do not connect the power cord immediately. The freezer needs to be placed in its intended location for at least 24 hours before connecting the power to make sure it will operate normally.

Alarm

DW-86L419/579/729/829

Alarm Type	Tricolor lamp panel	Buzzer	A prompt box pops up.
High-low temperature alarm	Red	Y	Y
Power-off alarm	Red	Y	Y
Door opening alarm	Red	Y	Y
Main temperature sensor fault alarm	Yellow	Y	N
Heat exchanger fault alarm	Yellow	Y	N
Condenser dirty alarm	Yellow	Y	N
High ambient temperature alarm	Yellow	Y	N
Battery fault alarm	Yellow	Y	N
Backup system fault alarm	Yellow	Y	N

Automatic recovery of alarm

• Under alarm, the beeper alarm could be ceased by pressing any key, and the fl ashing alarm and

remote alarm will not be ceased(DW-86L419/579/729/829:Click "×"on the upper right-hand corner in the prompt box to stop the buzzer alarm. If there is no prompt box,click the alarm on the bottom of the homepage to view the alarm, and then stop the buzzer alarm.).

• If the alarm condition is still available, the beeper alarm will be recovered automatically 30 minutes after ceasing.

Remote alarm terminal


• The remote alarm terminal is installed on the right side of the freezer's engine cabin, and the alarm signal will be output by this terminal. The load capacity of the terminal is DC 30V, 2A.

•Contact output:

The remote alarm terminal is divided into: Normally open, normally closed, common terminal.

Users may select normal open or normal close according to their own requirements.


 Standard R485 port can be used to transmit freezer temperature data to customer's software for monitoring

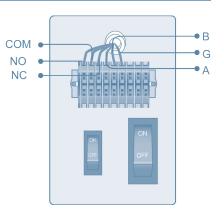

(î:	Help	2017-03-01 11:20:56 Administrator
	Inbox	Sent box
	title	Message sender:User1 2018-10-12 10:32
	Start time -	End time Inquiry Delete Select all

Notebook

It can be used to create new message, and view the records. For previous records, you can directly click them for modification and saving in the pop-up interface; when clicking [Delete], you can delete one or all messages.

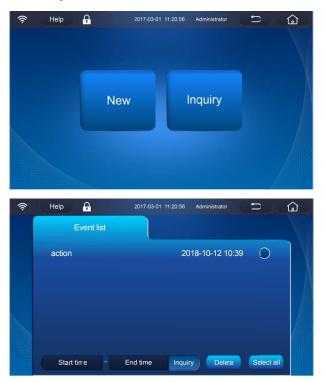

Initially Powering Up

When the unit is started for the first time, please follow the procedures below.

① While keeping the unit empty, plug in the power cord to a dedicated power socket that meets all requirements.

② Please connect the freezer to the power supply, turn on the protector located on the right of the freezer (as in the illustration on the right), and then turn on the battery switch. Protector Battery switch

③ If the unit has a backup cooling system (optional), turn off the backup system.

④ Set the unit to a desired temperature: Do not load the unit with any samples. Power up the unit to let it run down to -60° . Let it run at -60° for 8 hours then lower it to -80° . Observe the unit performance for 24 hours for normal cycling to ensure that it is working properly.

(5) Once the unit is confirmed to be operating properly, it is ready to be loaded with samples. In principle, the freezer unit should be set at about 3°C above the desired temperature. For example, if the storage temperature is -60°C, set the unit at -57°C. Load the unit with samples in batches of less than 1/3 of the unit's capacity. Make sure that the unit is capable of cycling for more than 8 hours.

(6) If the unit has a backup cooling system (optional), turn it on.

If the freezer temperature warms up, and the issue cann't be resolved within a short time, please transfer the stored products to another freezer to avoid potential damage..

• Before placing samles in the freezer for storage, check the freezer temperature to make sure the temperature meets the storage requirement for the samples. This is to prevent any potential damage to the samples due to incorrect storage temperature.

• Due to the fact that it would take time to remove the heat from the stored product, there can be a difference between the set temperature and the displayed temperature. This difference will get smaller as the set point is lowered.

 Ultra low temperature freezers are designed to store samples. Therefore, they are not suitable for heavy loading with a large amount of products at one time as the operation can cause premature compressors over heating. Samples should be loaded in small batches. The temperature setting should be lowered incrementally, allowing the unit and products are stablized at each setting until the final temperature is reached.

• Do not use any unauthorized mechanical tools or other means to accelerate the defrosting process.

- Do not tamper with the refrigeration circuit.
- Do not use any non manufacturer-approved electrical components in the freezer.

Operation after a Power Outage


Haier ULT freezers' control settings are stored in their memory system. Should there be a power outage and recovery, freezer units can resume its operation based on the previous settings.



• In the event of a power outage and recovery, be sure to wait for at least 5 minutes before turning the unit on again to avoid damaging the compressors and refrigeration system.

• If the unit is not in use for a long period of time, make sure the power cord is unplugged. Deteriorated insulation of the power cord may lead to electric shock or fire.

• If the freezer is not in use in an area without any supervision, please make sure children will not approach the freezer and the doors should not be closed.

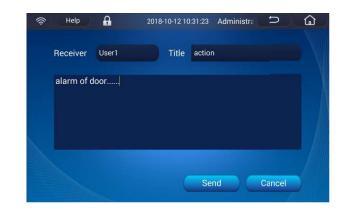

Message box

[Inbox]

You can click "Inbox" to view the messages that have been received. When a message has been read, its fonts become gray.

Inbox 💶	Sent box
title	Message sender:User1 2018-10-12 10:32
Start time -	End time Inquiry Delete Select all

[Sent box]

You can click "Sent box" to view the message that has been sent .

The user can query the name, model and machine code of this freezer, as well as version information of touch screen program.

You can view the electronic manual by clicking [Electronic manual].

Configuration: Click "Download" to download the set parameters of this freezer to a USB disk in the established format via USB, and then plug the USB disk that has saved the downloaded file to the USB interface of other freezers; after that, click "Upload" to upload this configuration file to other freezers, so as to realize that multiple machines can share the set parameters.

Database: Click "Download" to download the temperature data, alarm records, event records of this freezer to a USB disk in the established format via USB, and then plug the USB disk that has saved the downloaded file to the USB interface of other freezers; after that, click "Upload" to upload this database to other freezers, so as to realize that multiple machines can share the data.

Upload/download configuration files or databases: They can be uploaded or downloaded optionally via USB, so as to achieve the purpose that multiple machines can share the set parameters or saved data.



•When export data to USB disk, please use an USB disk with a file format of FAT32. •If the FAT32 USB disk cannot export data, please backup the file of the USB disk first and then format the USB disk or try another brand FAT32 USB disk.

5.Message




•New message

Create a message by clicking on "New message". In the interface of "New message", click "+" to select the users whom the message will be sent to, enter a title in the title bar, and enter the message content in the text box, then click "Send".

Freezer parts and control panels

Freezer parts

•DW-86L419/579/729/829

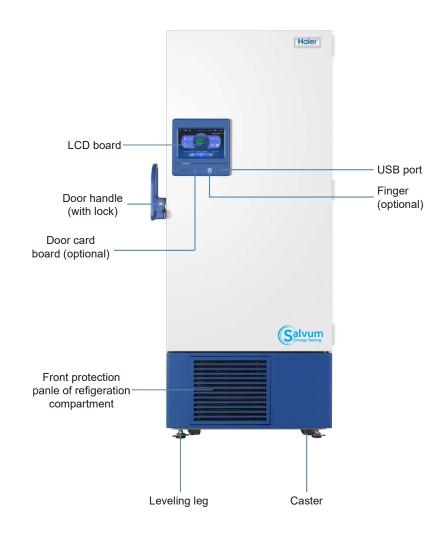

Control panel

•DW-86L419/579/729/829





For the accumulated injection time, it displays the accumulated injection time since the last start-up.



Enter/input "Start time-End time", and then click "Inquiry" to query the spraying record in this interval. One or several injection history can be deleted to the spraying history.

Device information

It can be used to query and save the related information, electronic manual of the freezer, as well as upload/ download its configuration file and database information.

📀 📀		2022-08-22 15:36:49 Login
	Name:	
	Model:	DW-86L829BPT
	Serial No.:	BE0F01EBA00QGN2P0005
	Version No.:	1.0.23.9.22082211_release
	Database:	Upload Export
	Setup file:	Upload Export
		Electronic manual
		27

🛜 Help 🔒	2017-03-01 11:20:56	Administrator	⊃்டு
-86 ℃			
Set point: -86°C			W
High temperature alar	m:-76°C		N N
Low temperature alarn	n : -96°C		
Ambient Temp : 21°C			
Voltage : 223V		w stage Condense mpressor fan 1	r Condenser fan 2
Temp. Recorder			
	Accumulated c	perating time (h):3	112

For the operation state of compressor / fan, "Green" indicates that it is running;" Grey "indicates that it is not running.

The electronic temperature recorder is used for collect the temperature inside the freezer alone. When clicking to enter [Temp. Recorder], you can view the recorded internal temperature, and when clicking [Download], you can download the temperature data visa USB.



· Backup system

It can be used to view the operating status, injection status, accumulated injection time, and injection history. It can set when connecting with backup system, but cannot set when there's no backup system and the display will show in gray.

Application method

DW-86L419/579/729/829

1.Homepage



In the home page, it can display the real-time internal temperature, ambient temperature, set point, input voltage, backup system state (optional), network state, electromagnetic lock state (optional), user login status, freezer health status, temperature graph, and message/ note book, etc.. •Icon of real-time internal temperature:

The icon reflects the health status of the freezer by changing colors:When the system is running normally, the color is "Green";When the system occurs a general alarm that does not affect the internal temperature, the color is "Yellow", for example, ambient temperature alarm;When the system has a serious alarm affecting the internal temperature, the color is "Red", for example, power failure alarm.

- Click to enter the sample management login interface
- Click to enter the graph viewing interface, where the temperature records can be downloaded
- **Q** Click to enter the information query interface
- Click to enter the freezer setting interface
- 중 Indicate that the network is enabled
- G Indicate that the electromagnetic lock has been installed, which is in the locked state
- **P** Click to enter the interface of configuring the sample bank server address and port.
- Click this button to enter the Alarm Record interface and the alarm record can be downloaded. When occurring a serious alarm that affects the temperature inside the freezer, the icon turns red, and when occurring a general alarm not affecting the internal temperature, the icon turns yellow. The buzz can be cancelled after clicking.



Click to enter the interface of message and note book .

Backup system
With this icon/identification:Indicate that the system is turned on;

Green light :Indicate that it is being sprayed;

Grey light: Indicate that there is no spraying.

System A 🔵 When the A/ B system is not running, the icon will display "Gray"; when in normal

System B • operation, the icon will display "Green";and when a certain system fails, the icon corresponding to this system will display "Red". (only for dual-system products) Help: vou can click this button to enter the help interface.

Log in/ Log out: Click the Login button to pop up the user login interface,, and click the Logout to log out the current account number.

2.Setting

The page has six buttons: Temperature, Initial, Operation, Backup, Network and User.



•Temperature

It can set the internal temperature and high- / low- temperature alarm value. Select the internal temperature value or high-/ low- temperature alarm value to be set, and then use the button of "+"/"-" for adjustment(You can also pop up a value selection box by clicking the value, and select the value by scrolling the selection box up and down). Click Save to finish settings.





Alarm record

It can be used to query the login alarm, door alarm, power alarm, battery alarm, sensor alarm, temperature alarm, condenser alarm, and backup alarm.

All records	Login	Door opening	Power	Battery	Sensor failur
2017-03-0	01 11:20:56		Conden	iser sensor1 fa	ault alarr
2017-03-0	1 11:20:56		Conden	iser sensor1 fa	ault alarr
2017-03-0	1 11:20:56		Back	kup system se	nsor fau
2017-03-0	01 11:20:56		Conden	iser sensor1 fa	ault alarr
2017-03-	01 11:20:56		Conder	nser sensor1 fa	ault alarr
2017-03-	01 11:20:56		Admir	nistrator Login	warning

Input "Start time-End time", and then click "Inquiry" to query the corresponding records within this interval; when clicking "Download", you can download the recorded events through USB.

Operating status

It can be used to query the freezer's real-time internal temperature , set temperature, high-/ low-temperature alarm value, ambient temperature, input voltage, cumulative operation time ; the compressor, fan, heating wire (dual-system model)'s working state; as well as the data recorded by the electronic temperature recorder.

When clicking somewhere on the graph, you can view the temperature value at a fixed point.

Input both the starting time and the ending time, and then click [Inquiry] to view the temperature graph in this interval; but if the interval is more than one month, the graph cannot be viewed.

Input both the starting time and the ending time, and then [Download] to download the temperature data within this interval through USB; but when you do not enter any time, all temperature data will be downloaded by default.

DATA can be exported in CSV or PDF, when you choose PDF, format at the interval between the start time and end time can not exceed 7days.

4.Inquiry

This page has six query buttons: Event record, Power and battery, Alarm record, Operating status, Backup system and Device information.



Event record

It can be used to query the records of door opening/closing, setting & modification, data upload/ download and account login/logout.

(íċ-	Help	A	2017-03-01 11:20:56	Administrator	Ð	ŵ
	All records	Door opening	Setting	Data	Login	>
	2017-03	-01 11:20:56		Administr	ator Login	
	2017-05-	01 11:20:56		Use	er1 Logout	
	2017-05-	01 11:20:56		U	ser1 Login	
	2017-03	-01 11:20:56		Administr	ator Login	
	2017-03	-01 11:20:56	Adm	iinistrator Create	a userU s e	
	2017-03	-01 11:20:56		Administra	ator Login	П
	Start t	ime – Enc	time Inquiry) 🦲	ownload	

Input "Start time-End time", and then click "Inquiry" to query the corresponding records within this interval; when clicking "Download", you can download the recorded events through USB.

· Power and battery

It can be used to view both the main battery and the backup system battery's input voltage, output voltage and remaining battery level

[Set point]

It is used to set the required temperature inside the freezer, with the adjustable range of -10 $^\circ\!C$ ~-86 $^\circ\!C$.

[Low temperature alarm]

It is used to set the alarm value of too low temperature inside the freezer, with the adjustable range of -15°C ~-99°C, which needs to be lower than or equal to the internal temperature value for 5°C. When the set value of internal temperature is changed, if the low temperature alarm value is not lower than or equal to the temperature value inside the freezer for more than 5°C, the low temperature alarm value will be automatically adjusted to "the set value of internal temperature - 5°C. If the low temperature alarm value is lower than or equal to the temperature value inside the freezer for more than 5°C, the low temperature alarm value is lower than or equal to the temperature value inside the freezer for more than 5°C.

[High temperature alarm]

It is used to set the alarm value of too high temperature inside the freezer, with the adjustable range of 0°C ~-81°C , which needs to be more than or equal to the internal temperature value for more than 5°C . When the set value of internal temperature is changed, if the high temperature alarm value is not more than or equal to the temperature value inside the freezer for more than 5°C , the high temperature alarm value will be automatically adjusted to "the set value of internal temperature +5°C . If the low temperature alarm value is lower than or equal to the temperature value inside the freezer for more than 5°C , the low temperature alarm value will not change.

Initial

It can set the language, date, time. Click any one value to pop up a selection box, and then you can select more values by sliding the values in the selection box up and down. When the setting is complete, click "Save", and then you can see that the time in the upper middle of the screen will be synchronized to the set value.



[Language selection]

It can set the language types (Chinese and English, selectable).

[Time adjustment]

Set the time value to the local actual time. The default value is read from the main control board. If there is no communication with the main control board, the last displayed time shall be the default value. If the time value disaccords with local actual time, it will affect the time display of the curve.

[Temperature scale]

Set the temperature standard, Celsius temperature standard $\,\,{}^\circ\!\!C$ and Fahrenheit temperature standard $\,\,{}^\circ\!\!F$.

Operation

It is used to set the alarm volume, screen brightness, Screen auto-lock, Auto-logout time, operational mode, account login mode, time delay to alarm, default display range of graph, time delay to start, ambient temperature alarm, time to door opening alarm, Update and reboot.

((r	Help 🔒	2017-0	3-01 11:20:56 Administrator		,
	alarm volume:	•			ц)»
	screen brightness:		•		-ờ́-
	Screen auto-lock:	5min	Auto-logout time:	5min	
	operational mode:	High performance	Energy saving Aut	horized	Ordinary
	Time delay to alarm:	30min	Default range for gr	aph:	2h
	Time delay to start:	1min	Ambient tempera	ture alarn	n: 35°C
	Time to door opening	g alarm:	1min Update	Rebo	ot
	Cance		(Save	

[Alarm volume]

The alarm volume can be adjusted by sliding it. On the left side: Minimum volume; On the right side: Maximum volume Click to enter "Mute" (I).

[Screen brightness]

The brightness of the liquid crystal screen can be adjusted by sliding it. On the left side: Minimum brightness; On the right side: Maximum brightness

[Screen auto-lock]

It can be used to set how long the screen will enter the sleep state for protection after having not clicked it, so as to extend the service life of the screen, with the default values of 5min,5min,10min, 15min, 20min, and never selectable.if you choose "never" the screen will enter the screen saver 5 minutes later. After the screen sleep, click anywhere on the LCD screen, it will automatically wake up and enter the homepage.

Choose "never" no operation within 5 mins enter screensaver.

[Auto-logout time]

It can be used to set how long the logged-in account will automatically log out after having not carried out any operation, so as to protect the account security, with the default values of 5min, 5min, 10min, 15 min and 20min (selectable).

[Operational mode]

High performance: It can provide high temperature uniformity and stability of the temperature inside the freezer, so as to store the sample with a very severe temperature control requirements on the internal temperature.

Energy saving: Compared with high-performance mode, it can save 10% of electricity, which will be used to store the sample with a not too severe temperature control requirements on the internal temperature.

When the electromagnetic lock is not used, the registered card can be used for account login on the login interface ;

Old Password:	Register Card
New Password:	
	CARD
	Please swipe the card on the card board.

[Delete]

Select an existing account, and then click it for login to delete the existing account.

3.Graph

It is convenient for viewing the graph of temperature inside the freezer, and it can be used to select the graph within the required time period and then download the temperature record. In addition, the graph that exceeds the alarm value of high- and low- temperature will be shown in striking red.



[2h] It can be used to quickly view the temperature graph of the last 2 hours, with the default value of 2 h(the default display range of graph can be set in the operation settings);

[24h] It can be used to quickly view the temperature graph of the last 24 hours

[7day] It can be used to quickly view the temperature graph of the last 7 days

[one month] It can be used to quickly view the temperature graph of the last 1 month

In the graph chart, the middle line is the line of the set temperature inside the freezer; the upper / lower line is high-/low- temperature alarm value respectively.

The graph can be enlarged by sliding the screen.



User: In authorization mode, normal user can only view, query and save the freezer's information, as well as use message and note book functions, but cannot set the freezer.

Administrator: In authorization mode, only with administrator permissions, can the settings be changed. With administrator permissions, you can not only use all the functions of the freezer, but also change the authorization mode into the normal mode.

Register fingerprint : When using a fingerprint lock (optional), click "Register fingerprint " to pop up the prompt for fingerprint registration, and then complete the registration of fingerprint on the freezer. After that, the user can open the door by using the fingerprint or log in the account in the authorization mode. When unlocking the lock, press the fingerprint directly; when logging in an account, select "Register fingerprint " on the login screen.

When the electromagnetic lock is not used, the registered fingerprint can be used for account login on the login interface under the authorization mode;



Register Card: When using a card-punching lock (optional), click "Register Card" to pop up the prompt for card registration, and then complete the registration of IC card on the freezer. After that, the user can open the door by punching the card or log in the account in the authorization mode. When unlocking the lock, punch the card directly in the unlocking zone; when logging in an account, select "Register Card" on the login screen.

[User login mode]

Ordinary: In this mode, any user can mostly use all the functions of the freezer, including setting (except user settings) and modifying some information, but can't use the message function. Under the normal mode, if a user account has been created, the mode immediately changes to the authorization mode as long as this account is logged in.

Authorized: The authorization mode can protect the data set, as only the administrator can access to change the settings. Other users can only access to check the temperature in the freezer and alarm information and use the notepad and message functions.

Mainly applicable to the freezer used by more than one person or when the requirement for safety management of the freezer is strict.

[Time delay to alarm]

It can set the time interval for alarming once again when the fault has not been eliminated after setting "Cancel alarm", including pop-up of alarm box and buzz ,with the default values of 30min, 30min, 60min and 120min(selectable).

[Default range for graph]

It can set the default time range of graph display on the homepage of graph, with the default values of 2h, 2h, 4h, 6h and 12h(selectable).

[Time delay to start]

It is used to set the start-up delay time of high temperature compressor after switching on the power supply, so as to prevent multiple refrigerators from starting up at the same time, otherwise the power may be larger, resulting in insufficient power supply. Its default value is 1min, with the adjustable range of 1~15min.

[Time to door opening alarm]

It can set the alarm time of open door, with the default value of 5min(1 ~ 30min, selectable).

[Update] When connect to wifi, click "update" to check if there is latest LCD screen program in the server background. If so, it will update automatically (just available for Chinese market so for).

[Reboot] click "Reboot" to Reboot LCD screen.

[Ambient temperature alarm]

It can set the alarm temperature value of ambient temperature , with the default value of 35 $^\circ\!\!C$ (30 $\,\sim$ 50 ,selectable).

Network

It can set the network opening or closing : After the network search is completed, click the corresponding network, and enter a password for login, then connect to the network successfully. It is recommended for the user to open and then connect the wireless network, so that the touch screen can be remotely upgraded. If there is no wireless network, you can open the hot spot of the mobile phone for network connection, so as to remotely upgrade the program of touch screen.

Network : On Off	Search	
Network name : Haier		
H60-L02	হ	
CU_7ysd		
CU_iTV_7ysd		
beifensxk		
FAST_07D8		
baso88	<u></u>	

Network name: connected successful wifi name.

CU_iTV_7ysd Signal strength Strong Security level WPA-PSK-CCMP+TKIP][WPA2-PSK Password:	-CCMP+TKIP][WPS][ESS]
Wifi password	
Display password	
Cancel	Connect
beifensxk	

Backup

It can set ON/OFF of backup system, type of backup system, and injection temperature, meanwhile can projected injection time. It can set when connecting with backup system, but cannot set when there's no backup system and the display will show in gray.

	2017-03-01 11:20:56 Administrator 🗂 🔓
Backup system:	on of
Backup system type:	LN2 CO2
Set injection temperature	e:
Projected injection time:	0 hour Injection test
Cancel	Save

[Backup system]

When the user uses a backup system (optional), click "On"; and then click "Off" after the completion of use.

[Backup system type]

Two types of liquid nitrogen(LN_2) and liquid carbon dioxide(LCO_2) are available for selection(by default of liquid nitrogen).

[Set injection temperature]

Set the temperature when the backup system starting spraying: For liquid nitrogen, the default temperature is -70°C, with the adjustable range of 0 ~-100 °C; and for carbon dioxide, the default temperature is -70 °C, with the adjustable range of 0 ~-70°C; meanwhile, the set spraying temperature shall not to be lower than the high-temperature alarm value. When the set spraying temperature is lower than the high-temperature alarm value, it can not be saved.

[Projected injection time]

When installing a backup system, the user can fill in the estimated usable time of the installed backup system. When need to confirm the remaining conditions after the backup system has been used, the user can query its cumulative spraying time, and then judge the usage conditions of the backup system through the difference between these two.

[Injection test]

When the temperature inside the freezer does not reach the set spraying temperature, you can click "Injection test" to simulate the spraying conditions, so as to confirm whether the spraying system can be normally used. If the displayed temperature inside the freezer≥ the set temperature, after pressing the button of "Injection test", the displayed temperature is unchanged, and the indicator light of backup system on the homepage is ON; at this time, control to open the solenoid valve for spraying, and stop the spraying 10 seconds later. If he displayed temperature inside the freezer < the set temperature, after pressing the button of "Injection test", the displayed temperature changes rapidly to the set temperature; at this point, control to open the solenoid valve for spraying, and stop the spraying 10 seconds later; after that, the displayed temperature will return back to the actual temperature inside the freezer.

User

It can add, edit, delete user accounts, as well as set the user account permissions.



